
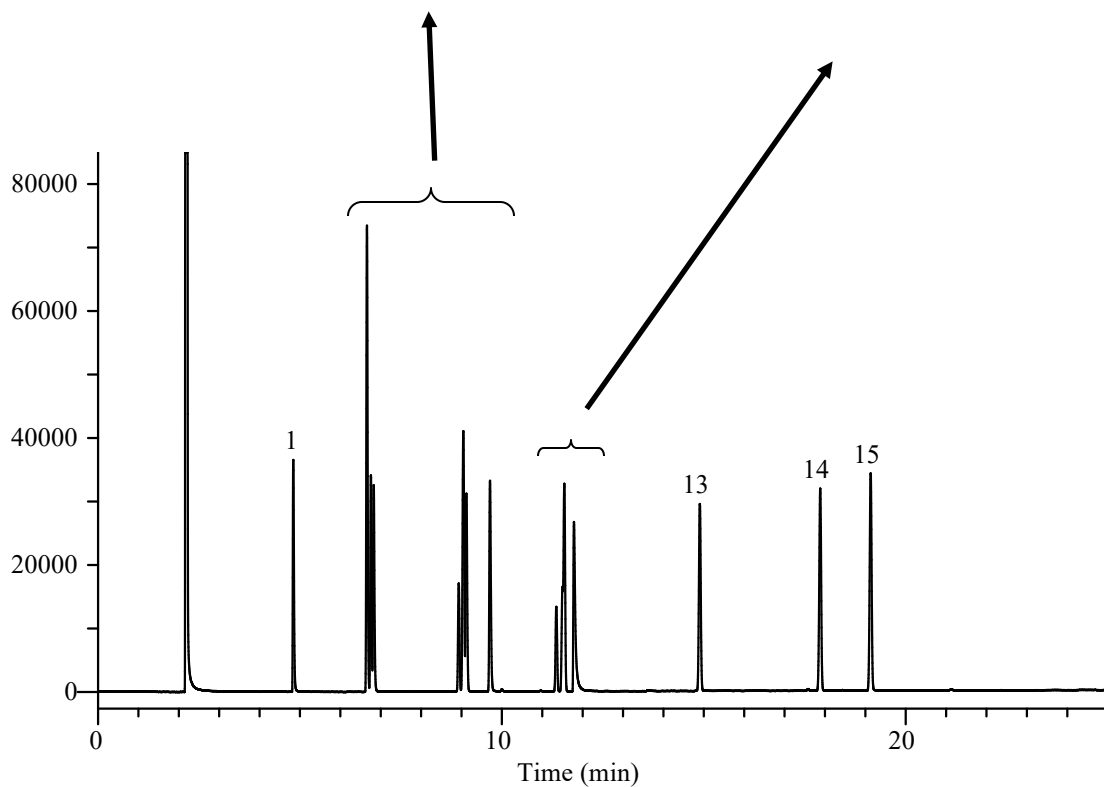


InertSearch™ for GC

InertCap® Applications

Flavor by Japanese Standards of Food Additives

Data No. GA173-0591

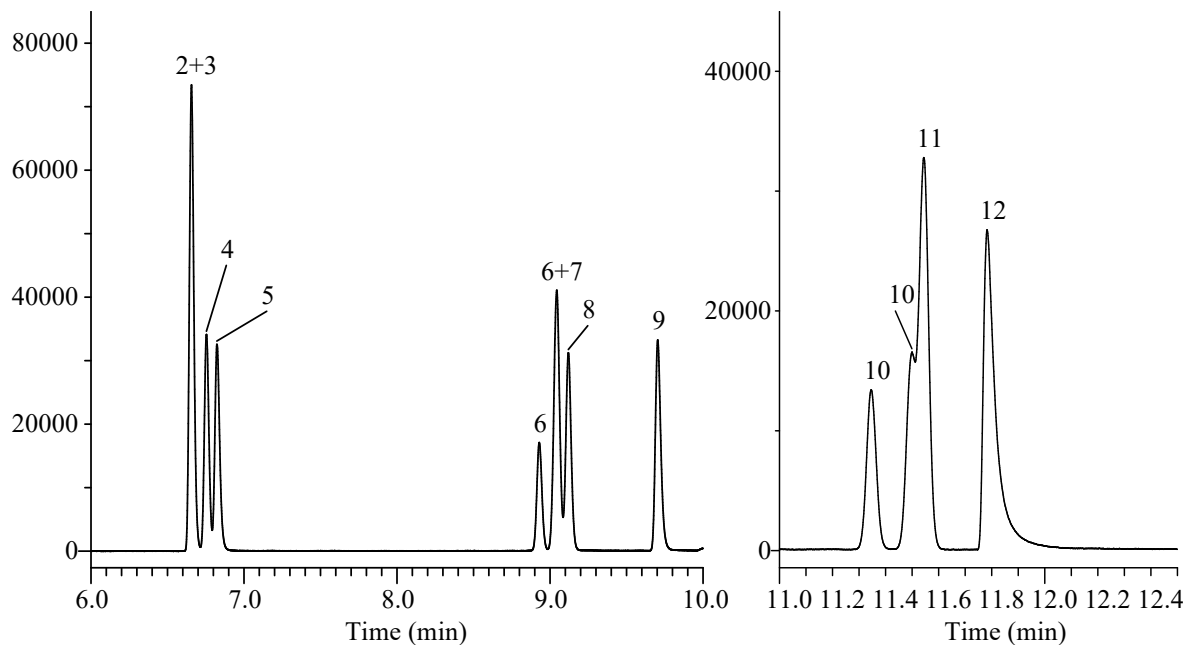

InertSearch™ for GC

InertCap® Applications

Flavor by Japanese Standards of Food Additives

Data No. GA173-0591

Conditions

System	: GC/FID
Column	: InertCap 1 0.25 mm I.D. x 30 m df = 0.25 µm
Col.Cat. No.	: 1010-11142
Col. Temp.	: 50 °C - 5 °C/min - 230 °C (4 min hold)
Carrier Gas	: He 80 kPa
Injection	: Split flow 115 mL/min 250 °C
Detection	: FID Range 10 ⁰ 250 °C
Sample Size	: 900 - 1100 µg/mL in Methanol 0.5 µL

Analyte

1. 2-Methylpyrazine
2. 2,5-Dimethylpyrazine
3. 2,6-Dimethylpyrazine
4. 2-Ethylpyrazine
5. 2,3-Dimethylpyrazine
6. 2-Ethyl-5(6)-methylpyrazine*
7. 2,3,5-Trimethylpyrazine
8. 2-Ethyl-3-methylpyrazine
9. 5-Ethyl-2-methylpyridine
10. 2-Ethyl-3,(5or6)-dimethylpyrazine*
11. 2,3,5,6-Tetramethylpyrazine
12. Phenethylamine
13. 5,6,7,8-Tetrahydroquinoxaline
14. 5-Methylquinoxaline
15. 6-Methylquinoline

*Isomers mixture