


Analysis of carbamate insecticides (Ditected by post-column method with OPA)

Data No. LA623-0626

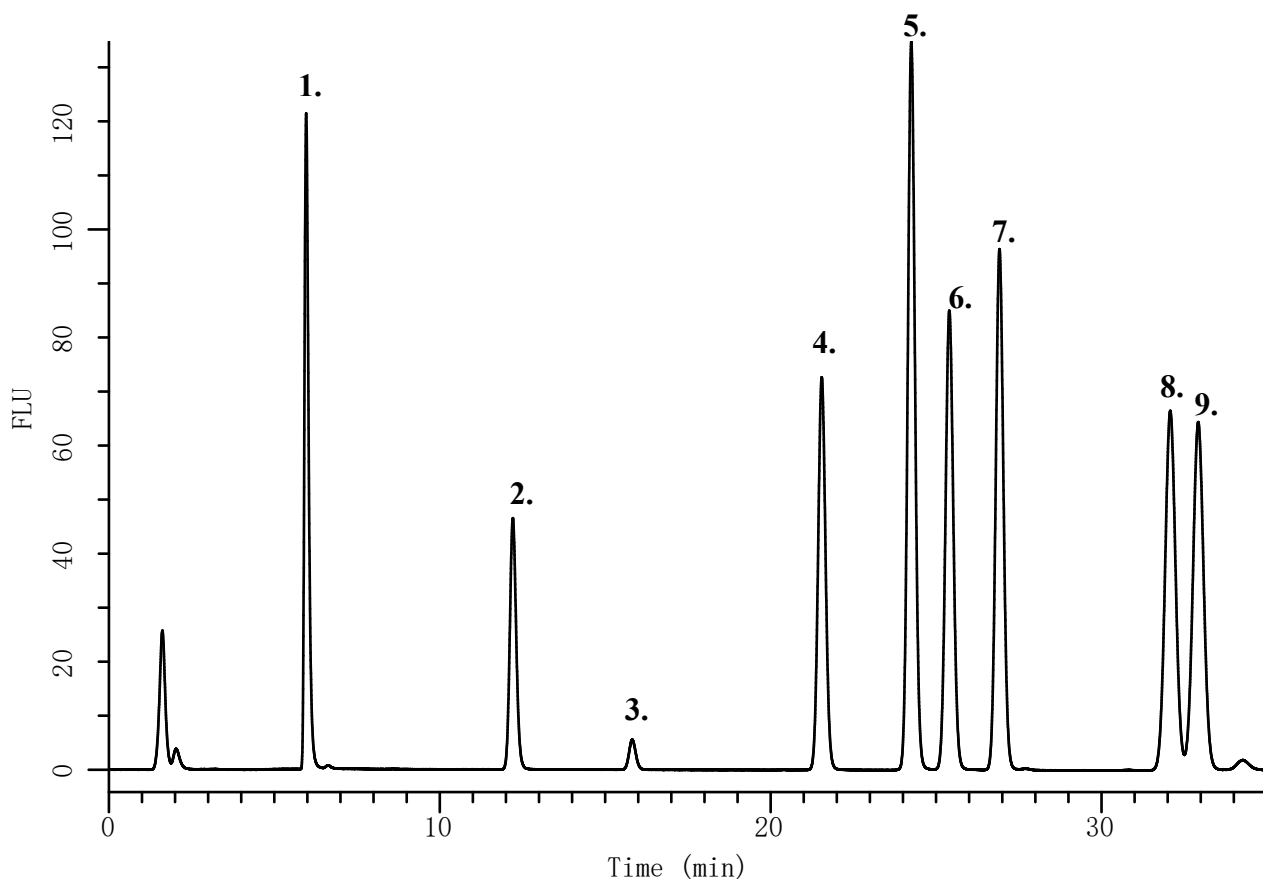

Conditions

System : GL-7400 HPLC system
Column : Inertsil ODS-3 (5μm, 150 x 4.0 mm I.D.)
Column Cat. No. : 5020-01761
Eluent : A) THF
 B) CH₃OH
 C) H₂O
 A/B/C = 0/12/88 - 0.1 min - 0/12/88 - 0.1 min - 10/0/90
 - 22.8 min - 33/0/67 - 12.0 min - 33/0/67 - 0.1 min - 0/12/88
 - 9.8 min - 0/12/88

Flow rate : 1.0 mL/min

Reaction Reagent : OPA発蛍光液*
 (Post-Column Derivatization)

Col. Temp. : 40 °C

Detection : Ex 339nm Em 445 nm (GL-7453A FL Detector)

Injection Vol. : 100 μL

Sample : Pesticides

* 発蛍光溶液: o-フタルアルデヒド10 mg及び2-メルカプトエタノール5 μLに
 0.05 mol/Lホウ酸ナトリウム溶液を加えて100 mLとしたもの。

Analyte:

1. Oxamyl	(1.0μg/mL)
2. Aldicarb	(1.0μg/mL)
3. Pirimicarb	(1.0μg/mL)
4. Bendiocarb	(1.0μg/mL)
5. Carbaryl	(1.0μg/mL)
6. Ethiofencarb	(1.0μg/mL)
7. Isoprocarb	(1.0μg/mL)
8. Fenobcarb	(1.0μg/mL)
9. Methiocarb	(1.0μg/mL)